Hinweis zur Montage der Membrane

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Notes regarding the installation of the pump diaphragm

Pump diaphragm/valve sets are designed to fit several models of diaphragm pumps, even though they have different performance specification.

Please follow all instructions regarding the replacement of diaphragms and valves listed in the operation manual. The service kit instructions may have changed since the operation manual was printed.

Compare the diaphragm with the one being replaced. If the original diaphragm has a square hole, simply punch out the section in the center of the diaphragm to convert it to a square hole and install as normal.

Attention: *Double diaphragm! Put the two diaphragms together with the printed sides outwards.*

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
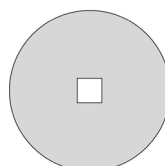
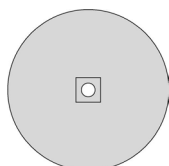
Attention: *Use the diaphragms only in the predetermined combination if the diaphragms are delivered packed in pairs.*

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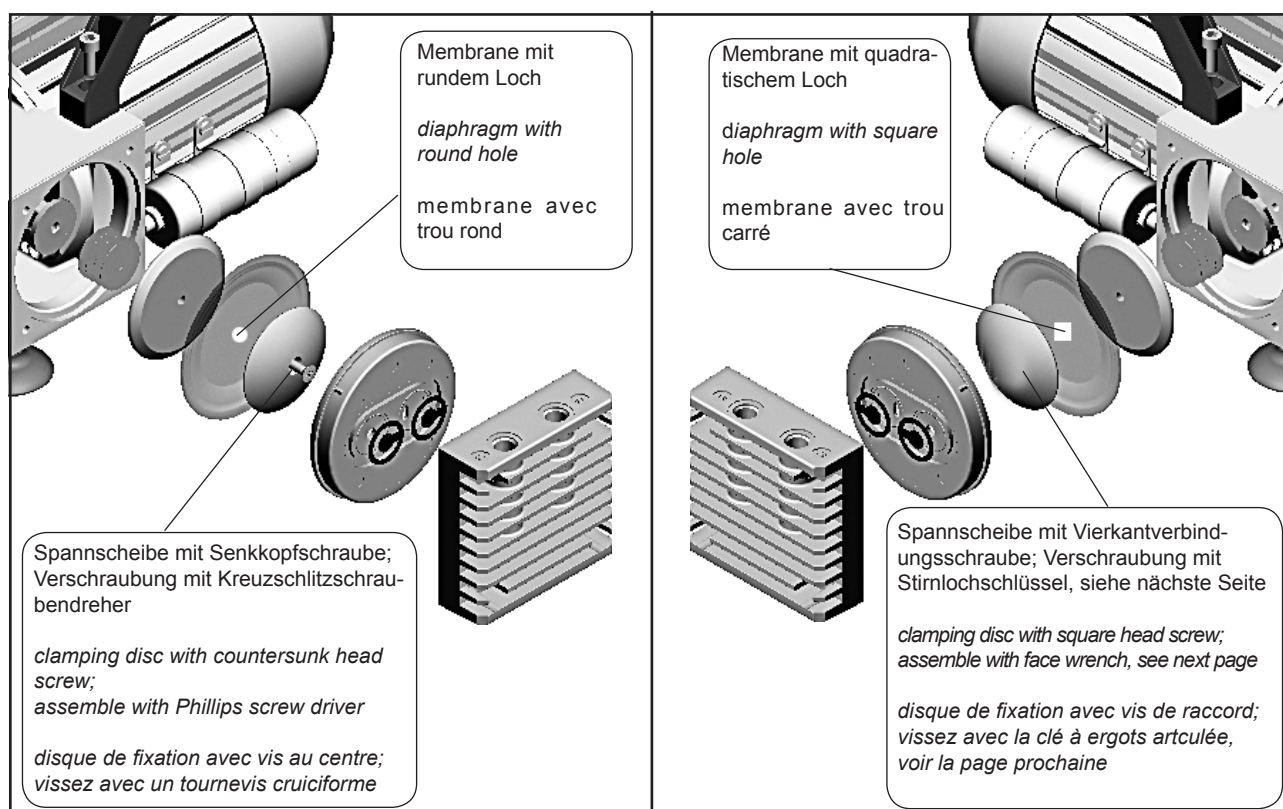Notes regarding the use of the "face" wrench

! Please follow all instructions regarding the replacement of diaphragms and valves listed in the operation manual. The service kit instructions may have changed since the operation manual was printed.

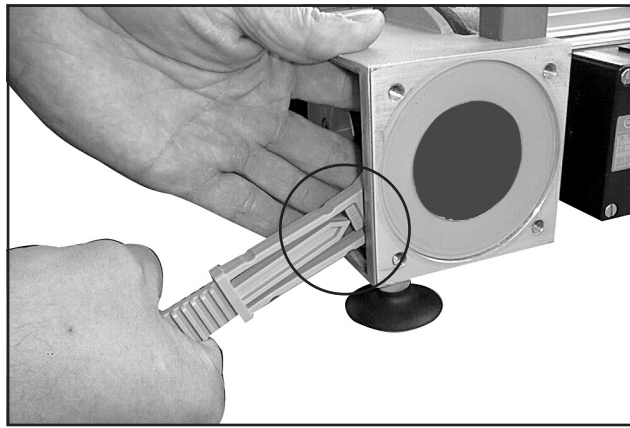
Use a face wrench with torque indicator (recommended: face wrench with torque indicator from VACUUBRAND, Cat.-No.: 20637580) to reassemble the clamping disc, diaphragm and tolerance washers and support disc to the connecting rod (see fig.).

Make sure that the square head screw of the diaphragm clamping disc is correctly seated in the guide hole of the diaphragm support disc.

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
Position 2: optimales Drehmoment

position 2: indicator should point here when the optimal torque is reached

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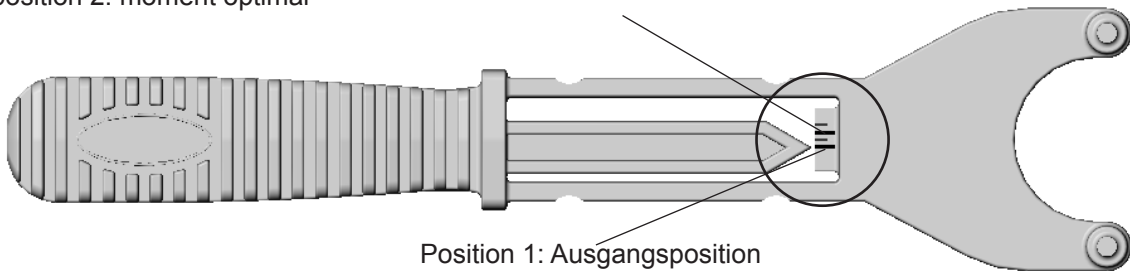*Optimum torque for the diaphragm support disc is **6 Nm**.*

The optimum torque is reached when the pointer in the handle of the face wrench points to position 2 which is the long line between the two shorter lines.

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Documents are only to be used and distributed completely and unchanged. It is strictly the users' responsibility to check carefully the validity of this document with respect to his product.

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