D

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GB

Notes regarding the replacing of diaphragms and valves

Comply with all instructions regarding the replacement of diaphragms and valves listed in the operation manual.

Before starting maintenance switch off the diaphragm pump and separate it from mains!

Attention: Double diaphragm (not MZ 2S NT)! Put the two diaphragms together with the printed sides outwards.

Attention: Use the diaphragms only in the predetermined combination if the diaphragms are delivered packed in pairs.

MZ 2S NT / ME 8S NT / ME 4R NT: Position diaphragm with pale side to diaphragm clamping disc (to pump chamber).

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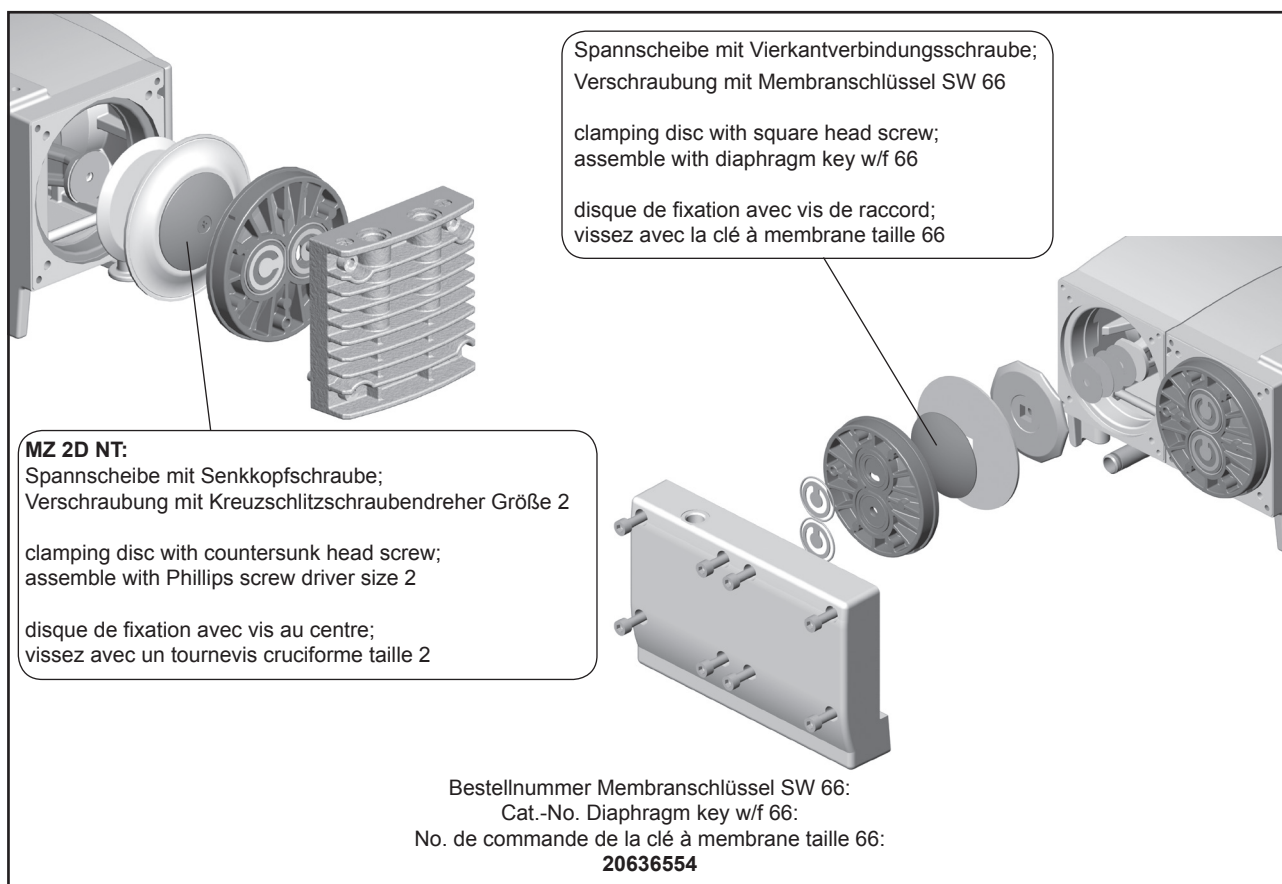
Documents are only to be used and distributed completely and unchanged. It is strictly the users' responsibility to check carefully the validity of this document with respect to his product.

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Assembling the diaphragm:

MZ 2D NT: Apply a drop of adhesive (OmniFit® 50M or Loctite® 243) to the lower side of the screw head and the thread and screw diaphragm clamping disc, diaphragm, diaphragm support disc and washers (if applicable) to connecting rod.

OmniFit® and Loctite® are registered trade marks of Henkel Technologies

Assembling with diaphragm key:

Position the new diaphragm between the diaphragm clamping disc with square head screw and the diaphragm support disc.

Make sure that the square head screw of the diaphragm clamping disc is correctly seated in the guide hole of the diaphragm support disc.

Lift the diaphragm at the side and position it carefully together with the diaphragm clamping disc and the diaphragm support disc in the diaphragm key (1).

Avoid damage of the diaphragm: Do not bend the diaphragm too much.

Screw diaphragm clamping disc, diaphragm, diaphragm support disc and washers (if applicable) to connecting rod (2).

Optimum torque for the diaphragm support disc: **6 Nm**, use torque key. Attach torque key to diaphragm key (hexagonal bolt 6 mm wide).

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© 2018 VACUUBRAND GMBH + CO KG Printed in Germany